

SPECIFICATION

DT-BLM-2038



DenimoTech

DT LABORATORY COLLIOD MILL

BENCH MODEL
FOR PREPERATION OF BITUMEN EMULSION
Sample size 500 - 3500ml.

Manufacturer:

*DENIMOTECH A/S
Haandvaerkervangen 12
DK-5792 Aarslev
Denmark*



PROCESS DESCRIPTION

The DENIMOTECH laboratory colloid mill – bench model, is especially designed for production of regular road grade bitumen emulsions that do not require production above boiling point for water.

The unit is one integrated unit consisting of the following parts:

- Colloid mill with specially designed high-speed AC motor.
- Thermostatically controlled electrical heating of colloid mill.
- Pipe system with funnel to receive the bitumen and other ingredients, re-circulation pipe system and sample valve.
- Electrical box with start/stop, thermostat, main power switch and internally; contactor motor protection and frequency controller.
- 4 litres funnel.

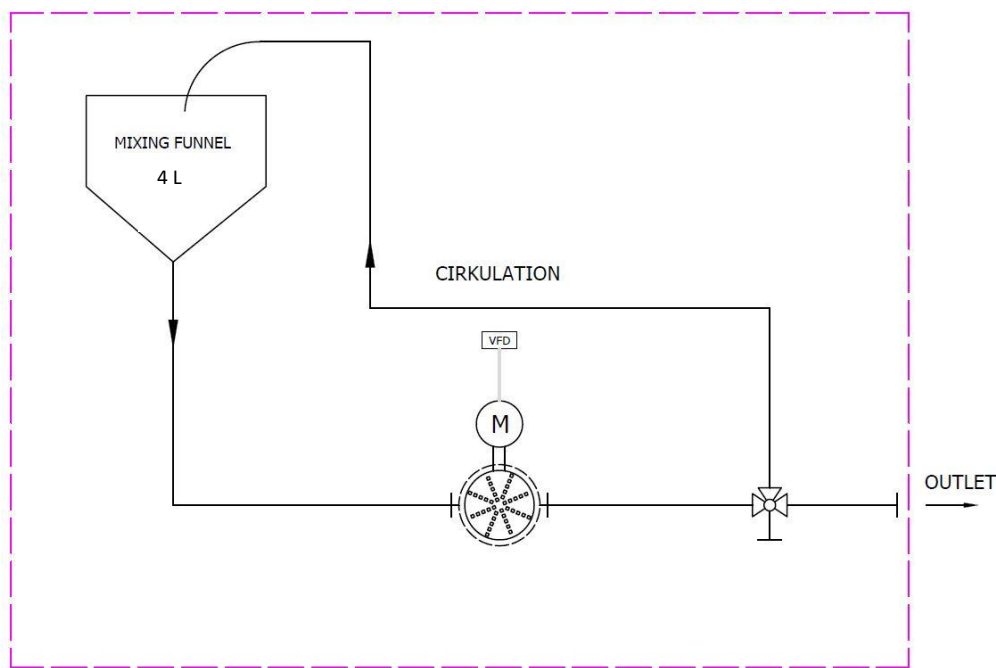


Figure 1: Ex. on a laboratory (bench model) plant flow-diagram / flow-process.

GENERAL FEATURES

Easy operation

The compact design of this unit including mill and motor, frequency controller and pipe system allows placement on a small frame that can be placed on a table and therefore easy to move around in the laboratory.

The electrical controls are collected in the control cabinet. From this you can operate the bench laboratory mill unit.

One person is enough to operate this unit.

Operational reliability

The plant is designed and build in full compliance with the latest European Standards of process machinery including mechanical and electrical design as well as documentation.

Only high-quality components and materials are used to ensure best possible operational reliability in tropical and tempered climates as well as industrial environment.

All parts are completely numbered and identifiable in manuals and part lists. Extensive manuals go with the plants and include process and electrical diagrams, maintenance and service instructions and technical data for the various components and systems. Repair readiness is achieved with a relatively small stock of spare parts.

Fast and easy installation

The DENIMOTECH bench laboratory mill unit are fully assembled with all the equipment. Calibration and testing are done at the factory prior to shipment.



DENIMOTECH colloid laboratory mill

The colloid mill is the same as in the DENIMOTECH Research plant and is of the same design as the DENIMOTECH full-scale production mills. The rotor and stator system are designed to enable achieving a fine and narrow particle distribution and the variable speed drive with a maximum of 11.500 rpm makes it possible to:

1. Vary and optimise the particle distribution and thereby the emulsion characteristics to adapt to prevailing requirements and to achieve close correlation with different full-scale production mills.
2. Using many different bitumen that normally form too coarse particle distributions resulting in poor emulsion quality.



All wetted parts of the colloid mill are made from AISI 316L and the pipe system and other components in AISI 316 stainless steel for corrosion resistance against low pH emulsions. The small volume of the pipe system minimises the raw material need and enables production of very small samples of 500 ml only for normally bitumen.

Production control panel

All operating buttons are placed on control cabinet.
The main switch is placed on the side of the cabinet.

Display for start/stopping the mill and speed control.

The signs along with the handles indicate which way to turn the valves.



TECHNICAL DATA

Plant model	DT-BLM (bitumen emulsion)
Sample size	500 – 3500 ml
Funnel capacity	4000 ml
Temperature, maximum	200°C
Viscosity	100 – 500 cSt
Motor power	2.0 kW
Motor speed, continuously variable	6.000 – 11.500 rpm
Rotor / stator gap, variable in steps	0.6 mm / 1.4 mm / 2.2 mm
Voltage	3 x 400V / 3 x 230V V, 50/60 Hz
Electrical data	
Total rated power	2,5 kW
L x W x H - approx.	500 x 300 x 650 mm
Weight - approx.	53 kg